

# Spiraltron™ Electron Multipliers

## Spiraltron™ Series of Electron Multipliers for High Pressure Applications



### Applications

- ✓ Elevated pressures
- ✓ Portable mass spectrometers
- ✓ Residual gas analyzers
- ✓ Vacuum loadlocks
- ✓ Small ID vacuum systems
- ✓ Array detections, such as magnetic sector mass spectrometers


### Features

- ✓ Longer life
- ✓ Higher dynamic range
- ✓ Reduced ion feedback
- ✓ High performance at elevated pressures
- ✓ Gains in excess of 100,000,000
- ✓ Noise less than 1 count/sec
- ✓ Excellent single ion sensitivity
- ✓ Small, compact linear configuration

PHOTONIS USA Spiraltron™ series of miniaturized channel electron multipliers can deliver high performance at pressures up to  $10^{-2}$  torr. The internal spiral structure facilitates low noise performance by significantly reducing the ion feedback typically generated from the high concentration of residual gas molecules in high pressure environments.

### Multichannel Construction

A Spiraltron™ detector is a Channeltron® electron multiplier consisting of six individual channels fed from a common collector cone. This multi-channel configuration provides a six-fold increase in surface area compared to single channel electron multipliers, leading to longer life. The spiral section of the detector consists of six single channels which are twisted – barber pole fashion – around a solid center. This geometry results in effectively six times the output surface area and, in addition, allows a straight channel geometry since the curvature preventing ion feedback is accomplished internally. This dramatically increases detector life and dynamic range.

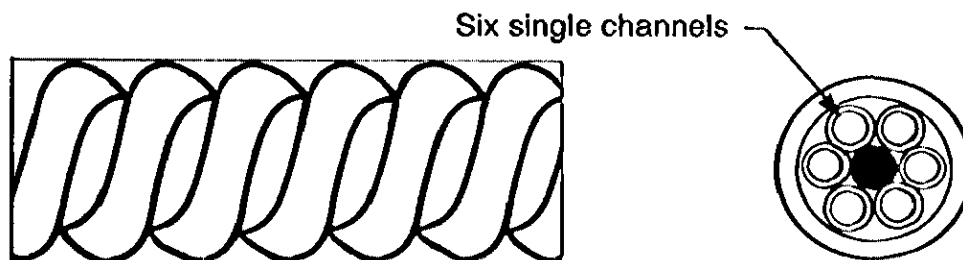

# Spiraltron™ Electron Multipliers

## Benefits of Spiral Channel Technology

- Compact, on-axis design - eliminates need for curved channels
- High dynamic range with six channels - delays the onset of saturation
- Higher pressure operation – spiral architecture significantly reduces ion feedback
- Analog or pulse counting modes
- Long Life - extracted charge comes from six channels, not just one.
- Cone shape can be tailored for the application - round, square, rectangular

## Spiraltron™

- Small, Linear Footprint
- Operates to  $1 \times 10^{-4}$  torr



Spiraltron™ detectors operate at elevated pressures into the  $10^{-4}$  torr range, a factor of 10 above CEMs and discrete dynode multipliers. Spiraltron™ electron multipliers are ideal for applications requiring high linearity. In addition, they provide excellent single ion sensitivity, producing Gaussian pulse height distributions (PHD).

## MegaSpiraltron™

- Compact and rugged design
- Operates to  $1 \times 10^{-2}$  torr



The MegaSpiraltron™ detector is a physically small, robust ion detector that can achieve high gain while maintaining low noise. These detectors have a compact, durable design and are only 1.35" long and 0.6" in diameter making them an excellent choice for portable instrumentation.

## MAGNUM™

- Easy integration
- Analog or pulse counting
- Operates to  $7 \times 10^{-4}$  torr



MAGNUM™ electron multipliers provide high dynamic range with high gain and low noise. The single piece cartridge design allows easy integration and replacement. MAGNUM™ detectors are available for analog or pulse counting applications, with standard or extended dynamic range performance.

*Channeltron, Spiraltron, MegaSpiraltron and MAGNUM are trademarks of PHOTONIS USA.*

The information furnished is believed to be accurate and reliable, but is not guaranteed and is subject to change without notice. No liability is assumed by PHOTONIS for its use. Performance data represents typical characteristics as individual product performance may vary. Customers should verify that they have the most current PHOTONIS product information before placing orders. No claims or warranties are made as to the application of PHOTONIS products. Pictures may not be considered as contractually binding. This document may not be reproduced, in whole or in part, without the prior written consent of PHOTONIS.