

Application note

Measurement of concentration of NH_3 in aqueous environment using integrated Rocket ATR spectrometer

Introduction

The Arcoptix OEM-ATR-20-12 spectrometer is an ATR (attenuated total reflection) version of the Arcoptix very successful range of spectrometers, adapted for measurements of liquids. This Application note is documenting the use of this spectrometer for measurement of diluted CO_2 in water.

Method

A 24% initial solution of NH_3 in water has been purchased. The solution was then sequentially diluted following a geometrical sequence with a factor of $\frac{1}{2}$. The final concentration in the sequence of measurements is in the range of 117 ppm.

Procedure:

- (1) The background spectrum was measured with degassed deionized water on the crystal, integrating individual 256 spectra with the resolution of 2 cm^{-1} .
- (2) The ATR crystal was then dried, and the sample was deposited using a micro-pipet.
- (3) The sample was left to stabilize for 1 minute and a spectrum integrating 256 spectra was taken in absorbance mode.
- (4) The recorded spectra were then investigated at 1101 cm^{-1} , where the maximum Absorbance of the diluted NH_3 is located.

Results

- (1) The detail of the absorption peak corresponding to the dissolved NH_3 for 1101 wavenumbers plotted for the variety of the concentrations is shown below.

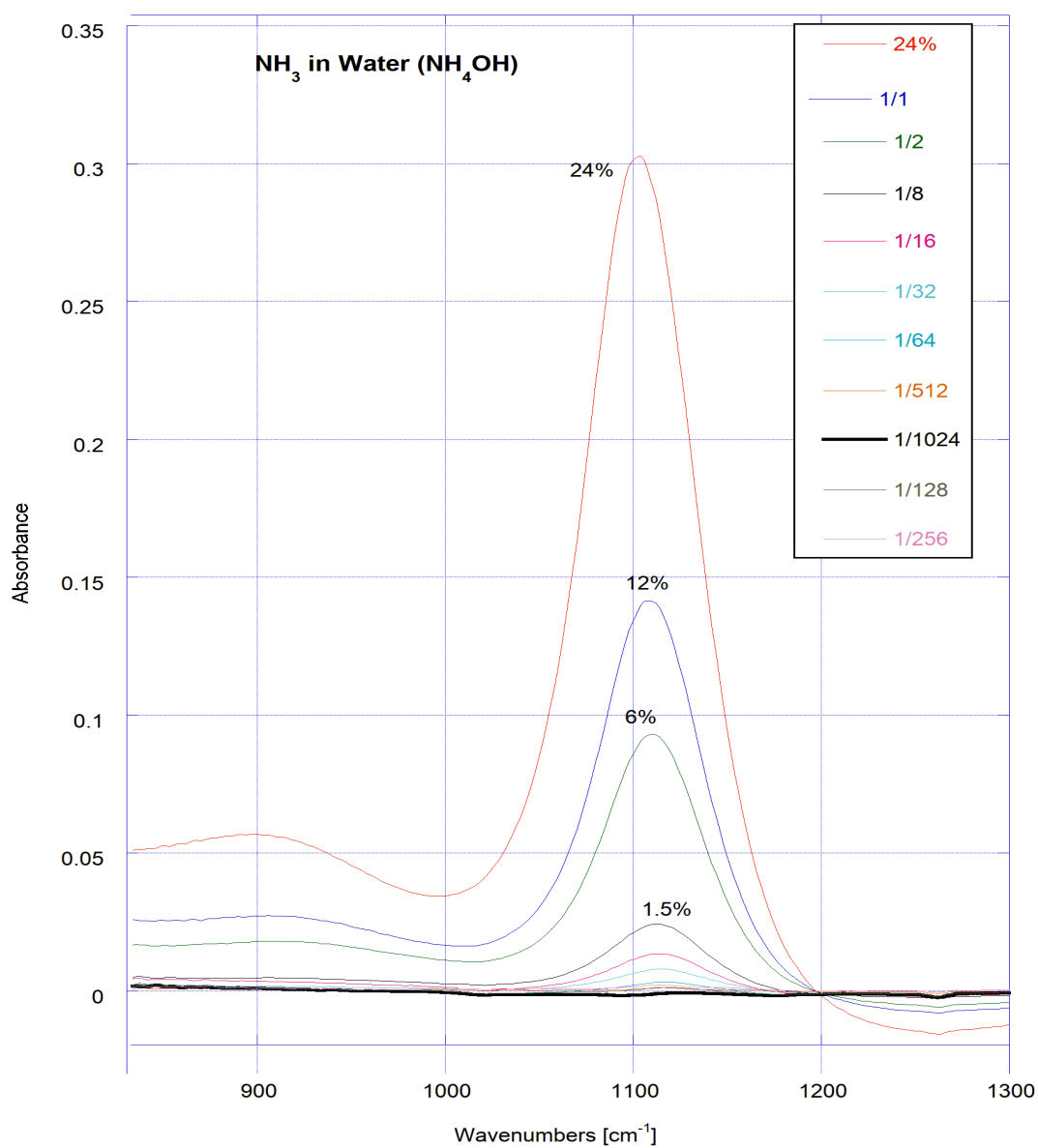


Figure 1 - CO₂ in water for various concentrations the dilution follows a geometric sequence with a factor of 0.5.

(2) A more detailed plot for the part of lower concentrations is shown below.

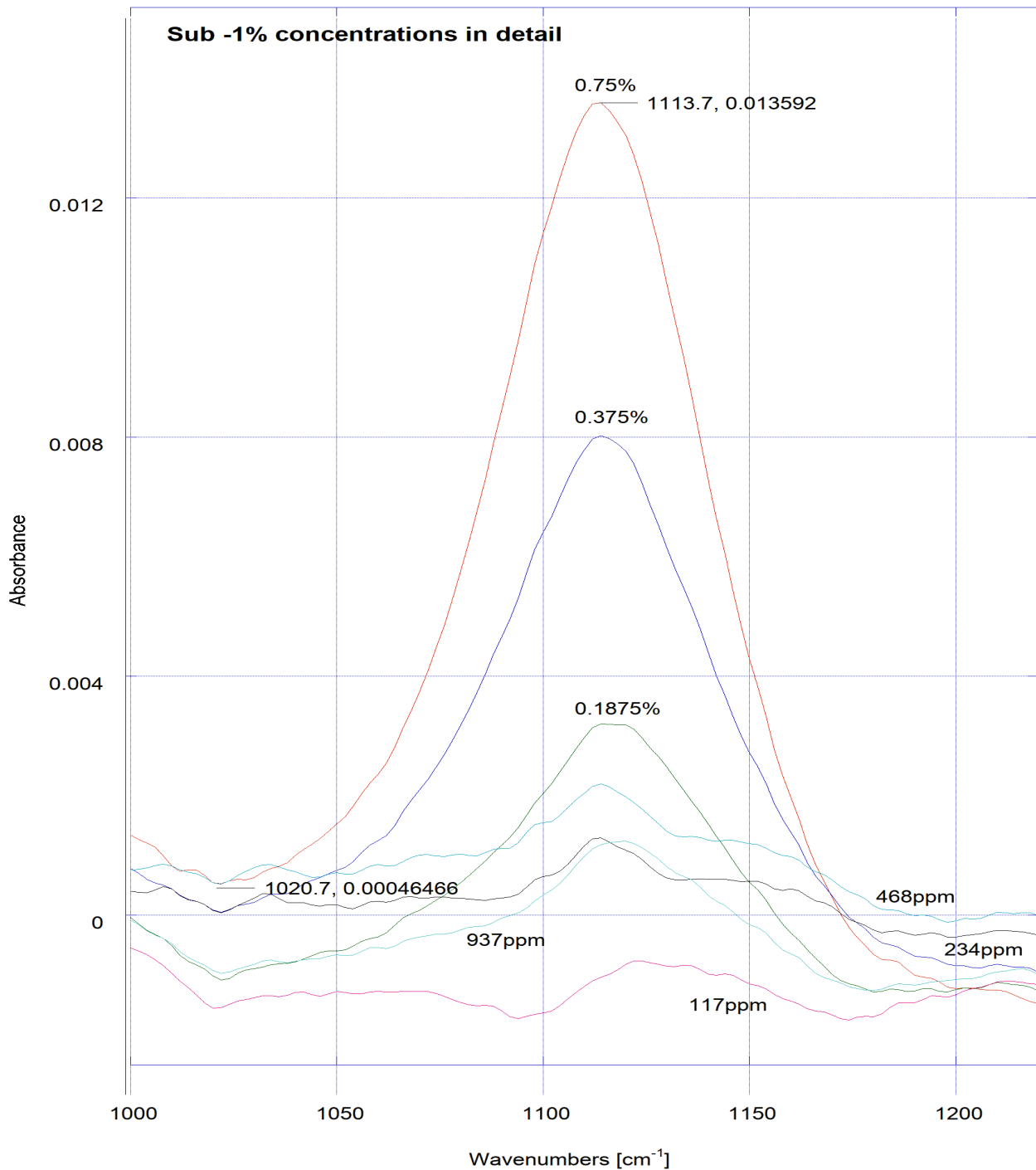


Figure 2 - Detailed plot for lower concentrations

(3) The sensitivity evaluation. For simplicity only the maximum of the absorbance signal with respect to the concentration at the 1101 wavenumbers have been plotted. With this respect the sensitivity of this method in the level of 1.1×10^{-6} of absorbance per ppm have been detected in the vicinity of the zero point (four lowest points have been line-fitted, and the slope has been evaluated). The LOD and LOQ determination would require more data in the dataset to be evaluated.

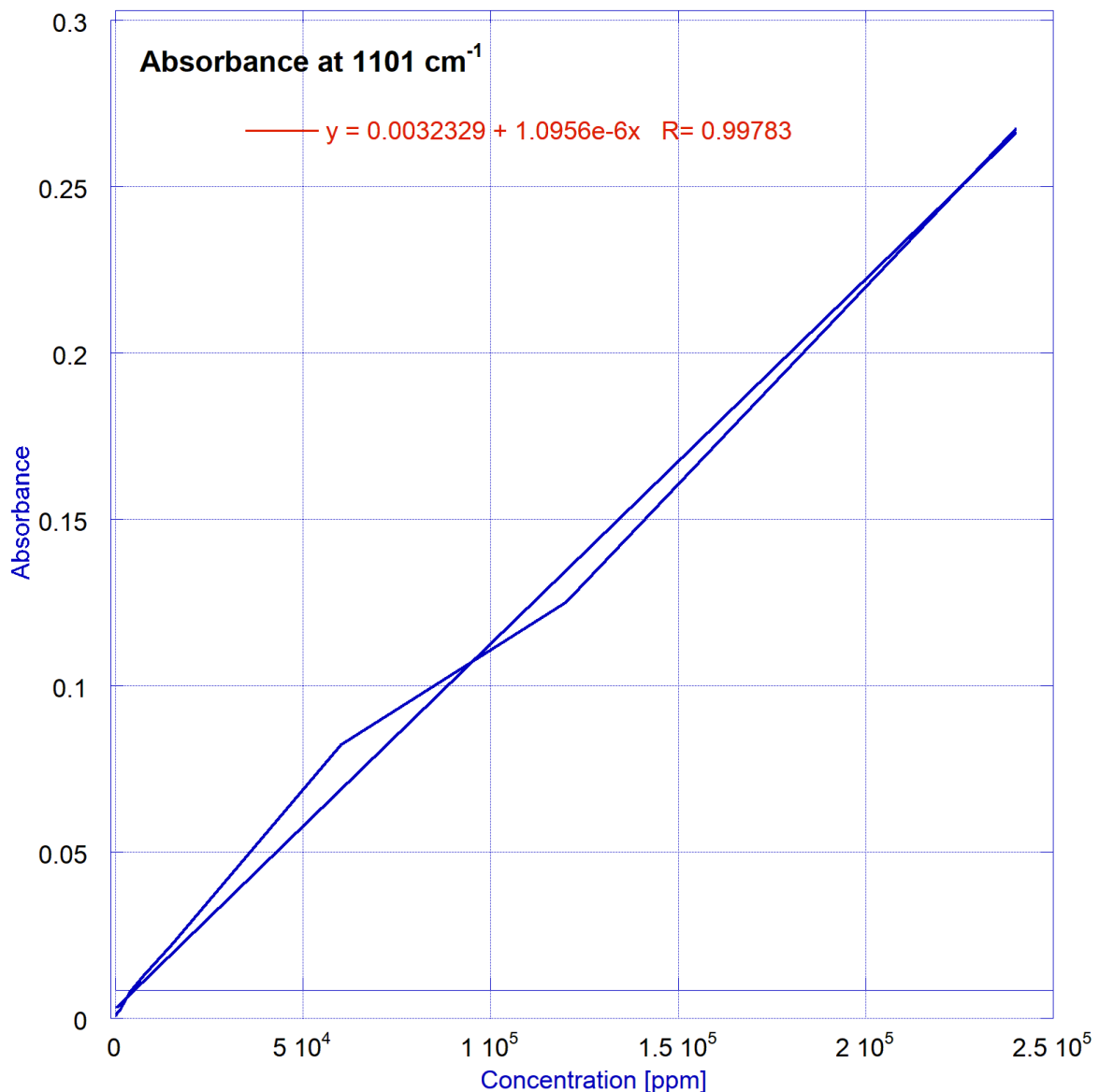


Figure 3- >Evaluation of the sensitivity of OEM- ATR-20-12 model for binary mixture of Ammoniac in water.

Arcoptix SA
Fabourg de la Gare 5A
CH-2000 Neuchâtel
Switzerland
Phone ++41 (0) 32 731 04 66
e-mail: info@arcoptix.com



Conclusion.

The OEM-ATR-20-12 has been evaluated as a suitable tool for measurements of the diluted Ammoniac in water on the 10 ppm level. Further – more sophisticated evaluation is to be performed to quantify the performance and to narrow down the margins of the precision.

 **東京インスツルメンツ**
TOKYO INSTRUMENTS
グローバルにネットワークを広げ、最先端の科学をお客様に提供

本 社：〒134-0088 東京都江戸川区西葛西6-18-14 T.I.ビル Tel.03-3686-4711
営業所：〒532-0003 大阪府大阪市淀川区宮原4-1-46 新大阪北ビル Tel.06-6393-7411
URL： <https://www.tokyoinst.co.jp> Mail： sales@tokyoinst.co.jp

 **Group Company**

 **UNISOKU**
TII Group

超高真空・極低温走査型プローブ顕微鏡
高速分光測定装置、クライオスタット

 **LOTIS TII**

Nd:YAGレーザー、Ti:Sレーザー
OPOLレーザー

- 本カタログに記載されている内容は、改良のため予告無く変更する場合があります。（製品の仕様、性能、価格などはカタログ発行当時のものです）
- 本カタログに記載されている内容の一部または全部を無断で転載することは禁止されております。
- 本カタログに記載されているメーカー名、製品名などは各社の商標または登録商標です。