

COMPACT VARIABLE LASER POWER ATTENUATORS – 990-0077, 990-0078

Features

- Attenuation range for transmitted beam:
 - Typical: 0.1% – 98%
 - For 257 nm and 266 nm: 1% – 96%
- Attenuation range for reflected beam
 - Typical: 2% – 99.9%
 - For 257 nm and 266 nm: 4% – 99%
- Extinction ratio: >1:1000
- Clear aperture:
 - 990-0077 – Ø17 mm
 - 990-0078 – Ø22 mm



990-0077 and 990-0078 Laser Power Attenuators

This manual variable attenuator series is designed to continuously attenuate linearly polarized laser beams. They feature precision opto-mechanics and high-LIDT optical components incorporated in a compact housing to ensure stable and reliable performance.

Attenuators are based on Brewster-type thin film polarizer fixed at 56° and a quartz half waveplate. The Polarizer's angle of incidence can be manually adjusted to get the best contrast or to manipulate the direction of the reflected s-polarized beam. The waveplate is housed in a 360° continuous rotation mount, which can be manually rotated to control the output power of the transmitted and reflected beams.


The intensity of either exit beam or their intensity ratio can be controlled over a wide dynamic range without alteration of other beam parameters. Pure p-polarization could be selected for maximum transmission, or pure s-polarization for maximum reflection. The s-polarization exit port can be blocked with a removable low-power dumper which absorbs laser power up to 5 W. External beam dumps are recommended for high-power applications.

990-0077 COMPACT VARIABLE ATTENUATORS WITH A CLEAR APERTURE OF Ø17 mm

Specifications

Clear aperture diameter	17 mm
Polarization contrast	>1:1000

Wavelength, nm	LIDT	Attenuation range (Tp)	Catalogue number
257	1 J/cm ² @ 257 nm, 10 Hz, 10 ns	1% – 96%	990-0077-257
266	1 J/cm ² @ 257 nm, 10 Hz, 10 ns	1% – 96%	990-0077-266
343	3 J/cm ² @ 343 nm, 10 Hz, 10 ns	0.1% – 98%	990-0077-343
355	3 J/cm ² @ 355 nm, 10 Hz, 10 ns	0.1% – 98%	990-0077-355
515	5 J/cm ² @ 515 nm, 10 Hz, 10 ns	0.1% – 98%	990-0077-515
532	5 J/cm ² @ 532 nm, 10 Hz, 10 ns	0.1% – 98%	990-0077-532
1030	10 J/cm ² @ 1030 nm, 10 Hz, 10 ns	0.1% – 98%	990-0077-1030
1064	10 J/cm ² @ 1064 nm, 10 Hz, 10 ns	0.1% – 98%	990-0077-1064

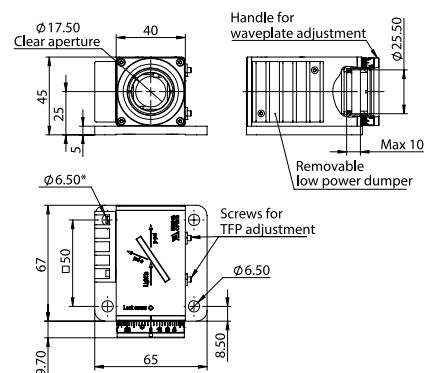


990-0078 COMPACT VARIABLE ATTENUATORS WITH A CLEAR APERTURE OF Ø22 mm

Specifications

Clear aperture diameter	22 mm
Polarization contrast	>1:1000

Wavelength, nm	LIDT	Attenuation range (Tp)	Catalogue number
257	1 J/cm ² @ 257 nm, 10 Hz, 10 ns	1% – 96%	990-0078-257
266	1 J/cm ² @ 257 nm, 10 Hz, 10 ns	1% – 96%	990-0078-266
343	3 J/cm ² @ 343 nm, 10 Hz, 10 ns	0.1% – 98%	990-0078-343
355	3 J/cm ² @ 355 nm, 10 Hz, 10 ns	0.1% – 98%	990-0078-355
515	5 J/cm ² @ 515 nm, 10 Hz, 10 ns	0.1% – 98%	990-0078-515
532	5 J/cm ² @ 532 nm, 10 Hz, 10 ns	0.1% – 98%	990-0078-532
1030	10 J/cm ² @ 1030 nm, 10 Hz, 10 ns	0.1% – 98%	990-0078-1030
1064	10 J/cm ² @ 1064 nm, 10 Hz, 10 ns	0.1% – 98%	990-0078-1064

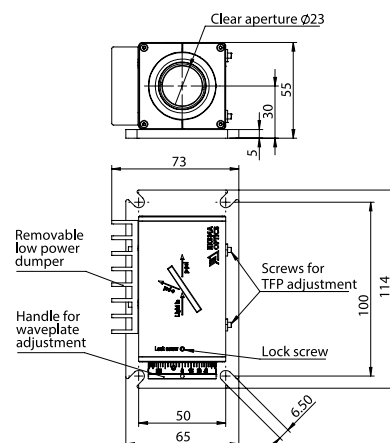



グローバルにネットワークを広げ、最先端の科学をお客様に提供

本 社：〒134-0088 東京都江戸川区西葛西6-18-14 T. ビル

Tel. 03-3686-4711

営業所：〒532-0003 大阪府大阪市淀川区宮原4-1-46 新大阪北ビル

Tel. 06-6393-7411

URL : <https://www.tokyoinst.co.jp> Mail : sales@tokyoinst.co.jp



超高真空・極低温走査型プローブ顕微鏡
高速分光測定装置、クライオスタット



Nd:YAGレーザー、Ti:Sレーザー
OPOLレーザー

- 本カタログに記載されている内容は、改良のため予告無く変更する場合があります。（製品の仕様、性能、価格などはカタログ発行当時のものです）
- 本カタログに記載されている内容の一部または全部を無断で転載することは禁止されております。
- 本カタログに記載されているメーカー名、製品名などは各社の商標または登録商標です。